

Technology that counts



NC-View™

Software User Guide

P/N 991-2022

Revision 1.12

30jan2023



ChemoMetec A/S

Gydevang 43 · DK-3450 Allerød · Denmark

Telephone: (+45) 48 13 10 20

Website: www.chemometec.com

E-mail: contact@chemometec.com

Caution!

This software must be operated as described in this user guide and documents referred to herein. Please read the entire guide and referred documents before attempting to use the software.

Contacting Support

Technical information that is not covered in this document or referred documents, is available from our support:

- Email questions to support@chemometec.com
- Speak to a Technical Support Specialist, by calling (+45) 48 13 10 20

Please create support files in NC-View™ via Help -> Create support files, before contacting ChemoMetec for support.

Sales and Ordering Information

To order NucleoCounter® NC-202™, NC-View™ software, and consumables, call (+45) 48 13 10 20 or send an e-mail to sales@chemometec.com.

Disclaimer Notices

The material in this document and referred documents is for information only and is subject to change without notice. While reasonable efforts have been made in preparation of these documents to ensure their accuracy, ChemoMetec A/S assumes no liability resulting from errors or omissions in these documents, or from the use of the information contained herein.

ChemoMetec A/S reserves the right to make changes in the product design without reservation and without notification to its users.

Copyright Notices

Copyright © ChemoMetec A/S 2020. All rights reserved. No part of this publication or referred documents may be reproduced, stored in a retrieval system, or transmitted in any form or by any means, electronic, mechanical, photocopying, recording or otherwise, without the prior written consent of ChemoMetec A/S, Gydevang 43, DK-3450 Allerød, Denmark.



ChemoMetec and NucleoCounter are registered trademarks owned by ChemoMetec A/S. Via2-Cassette and NC-View are trademarks of ChemoMetec A/S.


All other trademarks are the property of their respective owners.

Table of contents

QUICK GUIDE	4
GETTING STARTED	5
INTRODUCTION	5
GET STARTED IN EIGHT QUICK STEPS.....	5
STARTUP.....	6
VALIDATION WITH IQ/OQ PROTOCOLS	6
INTRODUCTION TO NC-VIEW™	6
NC-VIEW™ BASIC CONCEPTS AND TERMINOLOGY	6
USING NC-VIEW™	8
OPERATING THE MAIN WINDOW	8
USE THE CONTROL SECTION TO ACQUIRE DATA.....	8
<i>Sample and Dilution Volumes</i>	<i>9</i>
OPERATING THE IMAGE PANEL.....	10
<i>Configure Result Fields</i>	<i>10</i>
<i>Visual Inspection of Counted Cells.....</i>	<i>11</i>
FILE LIST IN MAIN WINDOW.....	12
<i>Tool Bar</i>	<i>12</i>
<i>Data in the File List.....</i>	<i>12</i>
<i>Define Columns to Display</i>	<i>13</i>
SELECTING A PROTOCOL	14
OPEN PREVIOUSLY ACQUIRED DATA FILES.....	15
EXPORT RESULT DATA	16
EXPORT PDF REPORTS	17
SOFTWARE OPTIONS	18
AUTO-EXPORT CSV FILES	18
AUTO-EXPORT PDF REPORTS	19
LICENSE(S) AND PROTOCOL LIMITATIONS	19
CERTIFICATE OF PRODUCT TESTING	19
INSTALLATION GUIDE	20
COMPUTER SPECIFICATIONS FOR RUNNING NC-VIEW™ SOFTWARE:.....	20
INSTALLING AND UPGRADING	20
UNINSTALL PROCEDURE.....	20
IMPORTING PROTOCOLS.....	21
MAINTENANCE AND BACKUP	21
TROUBLESHOOTING	22
CREATING SUPPORT FILES	24

Quick Guide

- ① Start the NC-View™ software via the *NC-View™* icon  in the Windows start menu.
- ② Once NC-View™ is open, NucleoCounter® NC-202™ will initialize. When the LED indicator on the instrument becomes green, the instrument is ready to use.
- ③ Select a protocol in the drop-down menu, or via the *Select Protocol* icon .
- ④ Load the Via2-Cassette™ with a cell sample and place it into the instrument cassette fixture.

OPTIONAL: enter *Sample ID* and *Operator name*
- ⑥ Start the analysis by selecting the *Run* icon  or press the RUN button on the instrument. A few moments later the result of the analysis will be displayed.

Repeat steps four, five and six to perform the same protocol on a new sample.


Getting Started

Introduction

The NC-View™ software is required for control and data acquisition with the NucleoCounter® NC-202™ instrument. NC-View™ is available for Windows 10 and 11 only – refer to the [Installation Guide](#) section for more information.



This manual is integrated into NC-View™. Pressing 'F1' while in NC-View™ will open the manual at the relevant section.

Get Started in eight quick steps

1. Log into your dedicated computer (see [requirements](#)) with your Windows administrator credentials.
2. Unpack the NucleoCounter® NC-202™ instrument and locate the software installation package on the USB key provided.
3. Launch the Install NC-View™ X.X.X.X.exe (the Xs indicate the version number e.g. 2.1.0.21.exe) using administrator permissions. **WARNING:** Do NOT open the .bin file.
4. Follow the on-screen instructions to install the camera and other device drivers. To complete installation, restart the computer.
5. Next, open the NC-View™  software in the Windows start menu.
6. Plug the included USB data cable into a USB 3 port on the computer.
7. Connect the other end of the USB data cable to the instrument using the screwdriver provided to secure the USB plug into the socket.
8. Insert the power cable into the outlet at the back of the NucleoCounter® NC-202™ and then connect it to a 3-prong outlet (i.e. including a grounding wire). The NucleoCounter® NC-202™ is ready to use when the RUN-indicator turns green (this may take about a minute).

For more details see the [Installation Guide](#) section below.




Startup

On the computer, open NC-View™ using the *NC-View™* icon  in the Windows start menu. The instrument will initialize, and the motor can be heard aligning during this process. When the LED indicator on the instrument turns green and the *Run* icon  becomes enabled, the instrument is ready to use. If the NucleoCounter® NC-202™ fails to connect, ensure that the instrument is connected to the computer using the supplied USB cable and ensure power supply is connected to the mains power outlet.

Validation with IQ/OQ Protocols

For the initial NucleoCounter® NC-202™ installation or when the instrument has been moved to a new location, performing a system validation is recommended using the *Installation Qualification* (IQ) and *Operation Qualification* (OQ) protocols.

Installation Qualification (IQ) and Operation Qualification (OQ)

1. Start NC-View™ using the *NC-View™* icon  in the Windows start menu.
2. Click the *Select Protocol* icon  to launch the [Protocol Browser](#).
3. In the Protocol Browser select the *NC-202 IQ/OQ* protocol.
4. Press the *Run* icon  and select *IQ* or *OQ* protocols, respectively.
5. Follow the on-screen instructions to perform the IQ and OQ test.
6. After the run has completed, a message will confirm whether the test has passed.

Introduction to NC-View™

For optimal user experience, a display resolution of 1920 × 1080 pixels or higher is recommended.

NC-View™ consists of multiple windows that can be moved and resized:

- [Main Window](#): Appears when NC-View™ is launched. It is used to operate the NucleoCounter® NC-202™, acquire data files and display acquired images.
- [Protocol Browser](#): Used to select and manage protocols.
- [File Browser](#): Shows the acquired data files (CM files) and is used to open files in the image display, export data or administer data files.

NC-View™ Basic Concepts and Terminology

- *CM file*: A ChemoMetec proprietary file format that contains the images, the result data and the metadata with file information. The file type is displayed as *.cm*
- *Protocol*: A file containing the settings on how the instrument acquires images, performs the image analysis and presents the results.


- *View*: The imaged area of a cassette. A view typically contains two images, one for the UV channel and one for the green channel. A CM file may contain a variable number of views depending on the protocol run.

Using NC-View™

Operating the Main Window



The Main Window is divided into three areas (indicated by numbers in the image below):

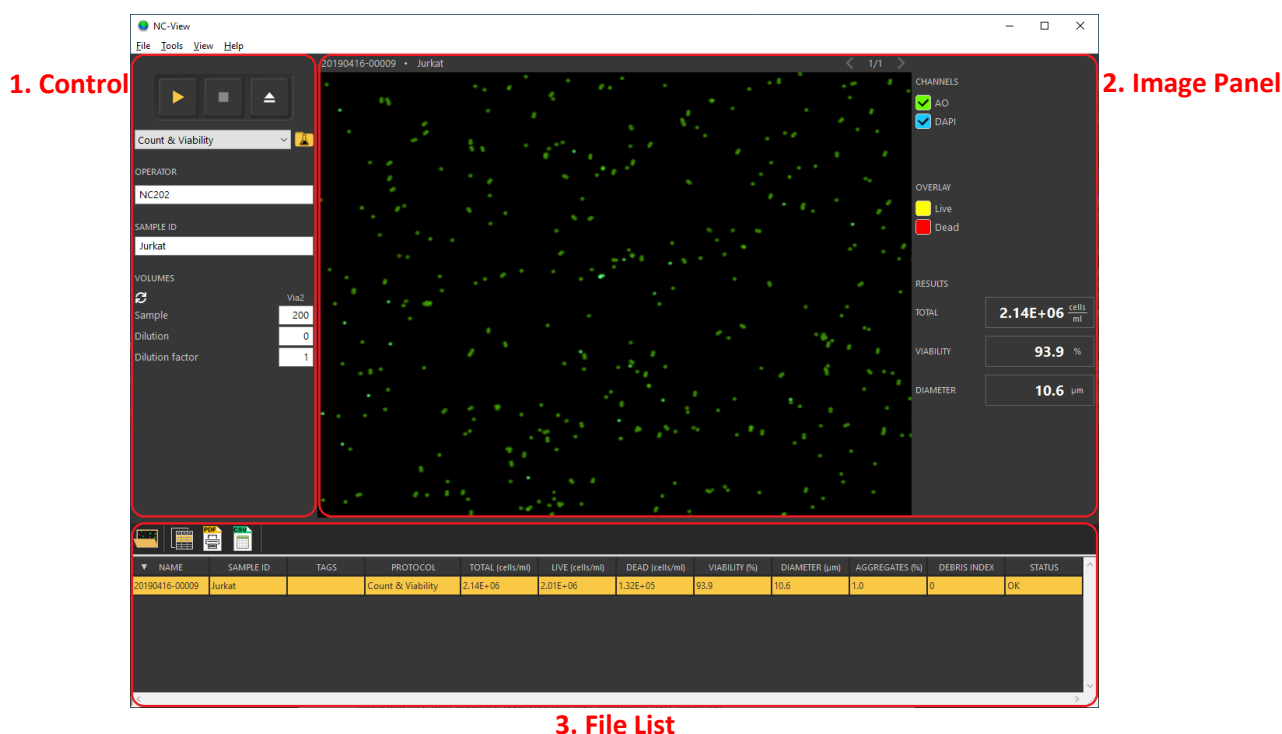1. The control section for [Acquisition](#) of new files.
2. The [Image Panel](#) displaying images and [Main Results](#).
3. The [File List](#) below the image populated with the files acquired on the same date as the CM file displayed.








Use the Control Section to Acquire Data

The control section on the [Main Window](#) is used for acquisition of new data.

1. Use the drop-down or the *Select Protocol* icon  and select the protocol of interest.
2. Load the Via2-Cassette™ with the cell sample and place it into the cassette fixture.
OPTIONAL: Enter *Sample ID* and *Operator name*.
3. Start the analysis using the *Run* icon  or the RUN button on the instrument.
4. The result will display after a few moments.


NOTE: Acquisition can be aborted using the *Stop* icon .

Once your analyses are complete, turn off the computer. There is no need to turn off the NucleoCounter® NC-202™ after use as the power consumption is very low when in standby mode. The instrument is calibrated by ChemoMetec and no subsequent calibration is required.

Graphical User Interface	Functions
	Run Start acquisition with the selected protocol.
	Stop Stop an acquisition session that is running.
	Eject Ejects the piston and releases the Via2-Cassette™.
Count & Viability 	Select Protocol Drop-down to select one of the ten last used protocols. Click on the folder icon to open the Protocol Browser , and select a protocol from the list.
OPERATOR NC202	Operator User-defined input: For multiple operators using the same Windows login.
SAMPLE ID <input type="text"/>	Sample ID User-defined sample naming field.
VOLUMES  Sample 200 Dilution 0 Dilution factor 1	Volumes The user can enter the volume before starting an analysis. The dilution or concentration of the sample will be factored in when calculating the results.


Sample and Dilution Volumes

Information about the sample dilution can be provided prior to starting an analysis, and the dilution or concentration of the sample will be factored in when calculating the results.

- When selecting a protocol, the volume fields will automatically update with the recommended volumes for the assay. The *Reset* icon  will reset the volumes to the default values.
- If the sample has been diluted, enter the sample and dilution volumes used, or simply enter the dilution factor.
- The dilution factor is defined as the sample volume divided by the sum of the sample and the dilution volumes. Entering a number greater than one in the input field represents a dilution of the sample, whereas entering a number less than one represents a concentration of the sample. Subsequently, the volume fields will automatically update to show the recommended volumes.
- If the sample has been concentrated, enter the original sample volume, and include the volume removed as a negative number in the dilution field or use the dilution factor as described above.
- Depending on the selected protocol, a solution field may be present. If the volume of the sample or dilution is changed and a solution input field is present, it will be updated with the recommended solution volume to add. If the user changes the volume in the solution input field, this volume will no longer be automatically updated.

Operating the Image Panel

Images of fluorescent cells can be viewed in the Image Panel. Use the mouse scroll wheel while the pointer is inside the image view area to zoom. To pan the viewed area when zoomed in, click and hold the left mouse button and drag to a new position. Display a CM file in the image display by clicking on the CM file in the [File List](#) or by opening the [File Browser](#) and double-clicking on a CM file.

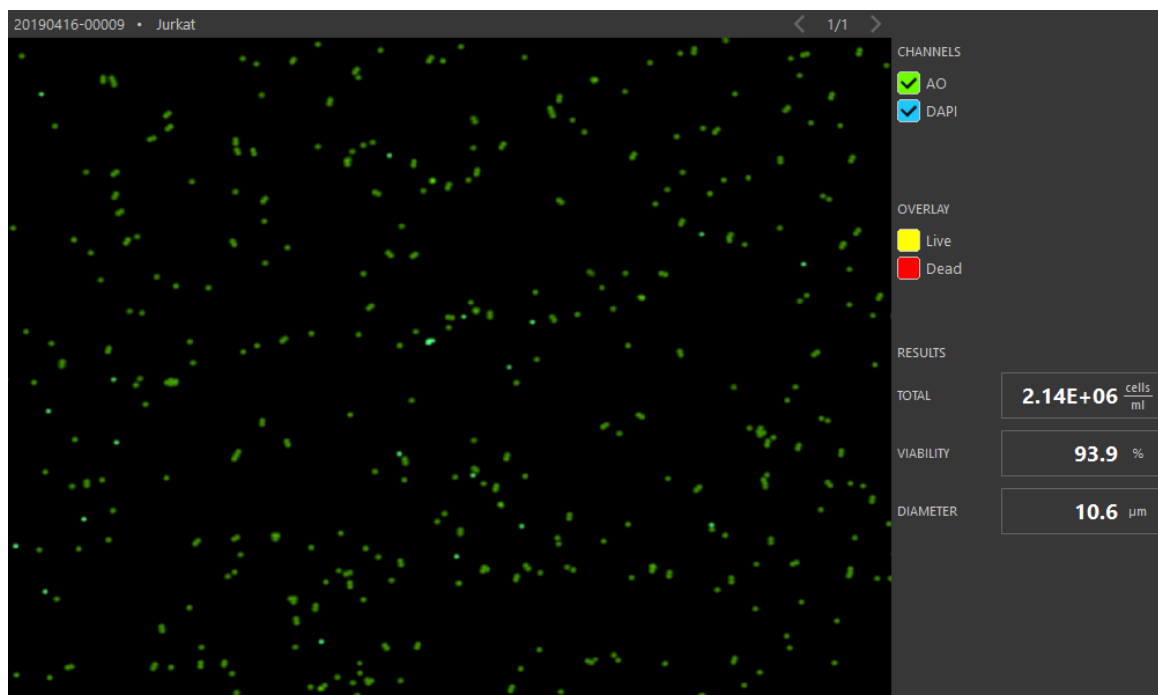


Configure Result Fields

On the right side of the image panel, three results fields are shown. These fields can be configured when a CM file is displayed by right-clicking on the results category, for example *Diameter*, and selecting the desired result output.

Graphical User Interface	Functions
20190416-00009 • Jurkat	Name and ID File name and sample ID of the current CM file displayed.
< 1/1 >	Previous and next view Click to change display if file contains multiple views.
CHANNELS <input checked="" type="checkbox"/> AO <input checked="" type="checkbox"/> DAPI	Channels List of channels in the current image. Click on a channel color to disable or enable the display of this channel.
OVERLAY <input checked="" type="checkbox"/> Live <input checked="" type="checkbox"/> Dead	Overlays List of overlays in the currently displayed image. Click on an overlay color to enable or disable the display of the overlay. Utilize this function for visual inspection of the counted cells.

Visual Inspection of Counted Cells

The overlays within a CM file can be used to visually inspect how the cells are counted by clicking the *Live* and/or *Dead* overlay checkboxes to enable the display of both or either overlays.

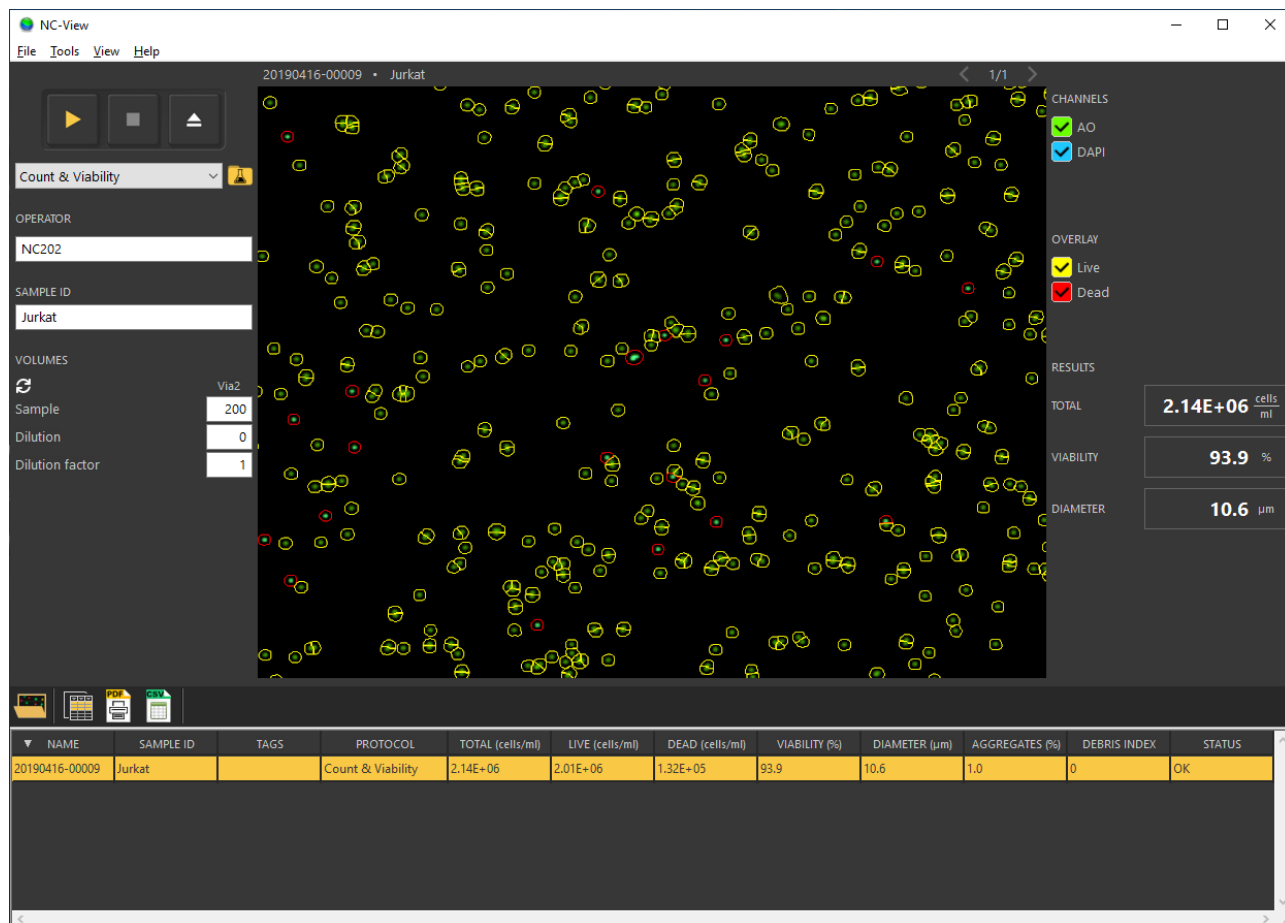

File List in Main Window

The file list below the image shows the files acquired on the same date as the displayed CM file.

Tool Bar

The tool bar at the top of the file list contains the following icons:



[File Browser](#): Find CM files that open in the image display or administrate data files.



[Export Result Data](#): Select file(s) and click the *Copy selected table data* icon to copy the data for the selected files to the clipboard. Data can subsequently be pasted into a spreadsheet.



[PDF Report](#): Create PDF reports for selected CM file(s).



[CSV Export](#): Create a CSV file for the selected CM file.

Data in the File List

There are two types of columns in the file list and [File Browser](#), *File information columns* and *Result columns*. The file information columns can be used to sort the file list by clicking on the column header. Sort the information in ascending or descending order by using the up and down arrows (▲ and ▼).

- File information columns
 - Name CM file name: date and run count.
 - Time Timepoint for when the CM file was created.
 - Sample ID User defined sample ID.
 - Tags Tags that have been added to the CM file.
 - Protocol Protocol name used when generating the CM file.
 - Media Media type (typically Via2-Cassette™).
 - User Windows username used when generating the CM file.
 - Operator Operator name used when generating the CM file.
 - Approved by Windows username used for approving the CM file.
 - Instrument Serial number of the instrument used for generating the CM file.
- Result columns
 - Total (cells/ml) Total cell concentration.
 - Live (cells/ml) Live cell concentration.
 - Dead (cells/ml) Dead cell concentration.
 - Viability (%) Percentage of live cells in the sample.
 - Diameter (µm) Median size estimation of the cells.
 - Aggregates (%) Percentage of cells in aggregates with five or more cells.

- Debris Index Index of debris events in relation to the cell count. For example, if the total cell count is 2×10^6 cells/ml and debris index is 50, the amount of debris is approximately 1×10^6 events/ml. The index value is rounded to the nearest 5.
- Status Shows *OK* or warnings such as *Foreign object detected* or *Out of range*


NOTE: Not all result columns will be available for every CM file type.

Warnings that appear in the status column for a CM file:

- *Out of range (5E+04– 1E+07 cells/ml)* The cell concentration in the Via2-Cassette™ is lower than 5×10^4 cells/ml or higher than 1×10^7 cells/ml
- *Bubble detected* A bubble that significantly affects the results is detected and removed from the analysis.
- *Foreign object detected* A foreign object that significantly affects the results is detected and removed from the analysis.



Define Columns to Display

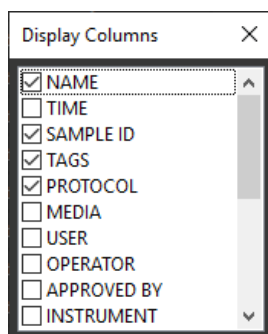
In the [File Browser](#) and file list in the [Main Window](#), columns can be shown or hidden by right-clicking on the heading row, opening the display columns dialog box and then selecting the column headings to be displayed.






NOTE: File *Name* is pre-selected


Selecting a Protocol

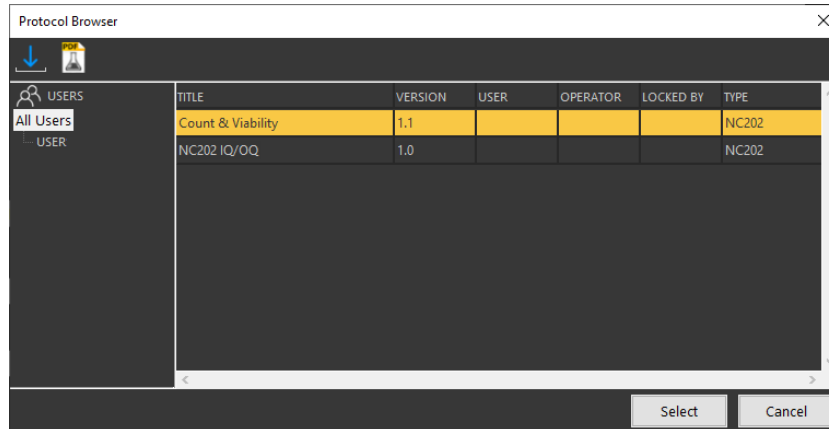
A protocol is a file that contains settings for how the instrument acquires images and how the images are analyzed. Protocols are selected via the Protocol Browser or the drop-down menu. The Protocol Browser is opened from the [Main Window](#) with the *Protocol Browser* icon . By selecting a protocol and clicking the *Application note*  icon, an application note for the selected assay will be displayed.



Graphical User Interface	Functions
	Application Note Opens the application note for the selected protocol.
	Import Protocol Opens a new dialog box in which cmsx or cmsu protocol files can be selected for import.
 USERS All Users USER	Users <i>All Users</i> or a single user can be selected as a search tool for locating the protocol of interest.





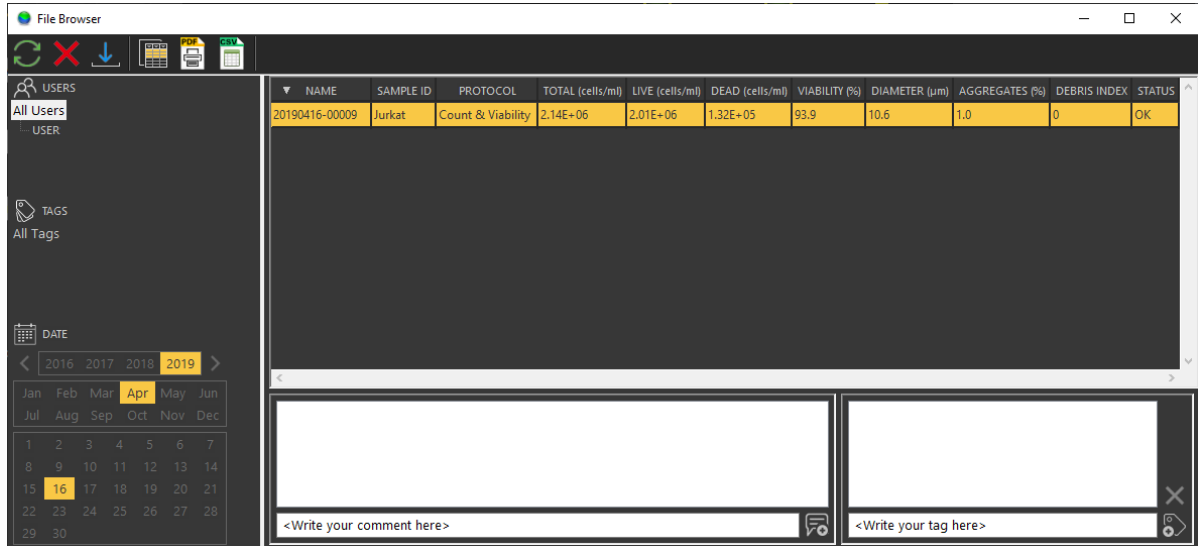







Open Previously Acquired Data Files

The File Browser provides an overview of the stored CM files. To open the File Browser, click on the *File Browser* icon  in the [Main Window](#). When double-clicking on a CM file, it will open in the image panel on the Main Window. This will also update the [File List](#) in the Main Window with the CM files acquired on the same date as the opened file. [Sorting](#) CM files, [Column Types](#) and [Column Display Configuration](#) work in the same way as with the [File List](#) (for more information see below).

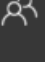


Right-click on one or more selected CM files to open a menu with the following options:



- *Reuse the run input parameters*: Updates the input fields on the Main Window with the parameters with which the selected file was acquired.
- *Copy file(s) to folder*: Opens a new window where files can be copied to another destination.
- *Sign Data*: When using the *Secure Mode* feature in NC-View™, a new window will appear, requesting authorization to sign the selected CM files.
- *Change sample ID*: The *Sample ID* for one or more selected files can be changed in a new window.
- *Properties*: File properties of a selected CM file open in a new window.

Icon	Functions
	Synchronize Synchronization will update the listing of the CM files in the File Browser.
	Delete file This action will delete the selected CM files.
	Import CM files Select the location of CM files to be imported into the File Browser. To export CM files simply select one or more files and drag and drop them to the desired destination.
	Copy selected table data Copies the selected table data to the clipboard so it can be pasted into another application.
	Create PDF report Select one or more files and click this icon to create a PDF report for each file.
	Create CSV file Select a file (one only) and click this icon to create a CSV file.
	Users Select all users or a single user to locate the specific CM files of interest.
	Tags Select all tags or a single tag to locate the specific CM files of interest.
	Calendar Select year, month, or day to locate the specific CM files of interest.
	Add Comment to CM file The user can add comments to a CM file. Note that comments are permanent and cannot be deleted.
	Add Tag User-defined keywords (e.g. project name) can be added to CM files as a tag for improved searchability.
	Delete Tag Select a tag in a CM file to delete.

Export Result Data


1. Select one or multiple files in the [File List](#) of the [Main Window](#) or in the [File Browser](#).
2. Click on the *Copy selected table data to clipboard* icon  to copy the data for the selected files to the clipboard.
3. Paste the data into a compatible application.

Export PDF Reports

1. Select one or multiple files in the file list of the [Main Window](#) or in the [File Browser](#).
2. Click the PDF report icon  to open the print options window.
3. The PDF report name is by default constructed as "<Instrument serial number>-<CM file name>.pdf". If a single file is selected, the PDF file name can be edited manually.
4. To select the destination folder for the PDF file(s), use the folder icon .
5. Select the print options (Merge PDFs, Add Sample ID to file name, Paper Size and format, Result Tables, Images, Signatures) and click *Ok* to save the PDF report(s).

NOTE: If Add Sample ID to file name box is checked, the sample ID will be added as a prefix to the PDF report name. The following characters are not allowed when this box is checked: * ? \ / " : < > | Each illegal character will be replaced by an exclamation point.

When selecting Merge PDFs for multiple files, the PDF reports are combined into one PDF document. The Add Sample ID to file name is not available with the Merge PDFs.

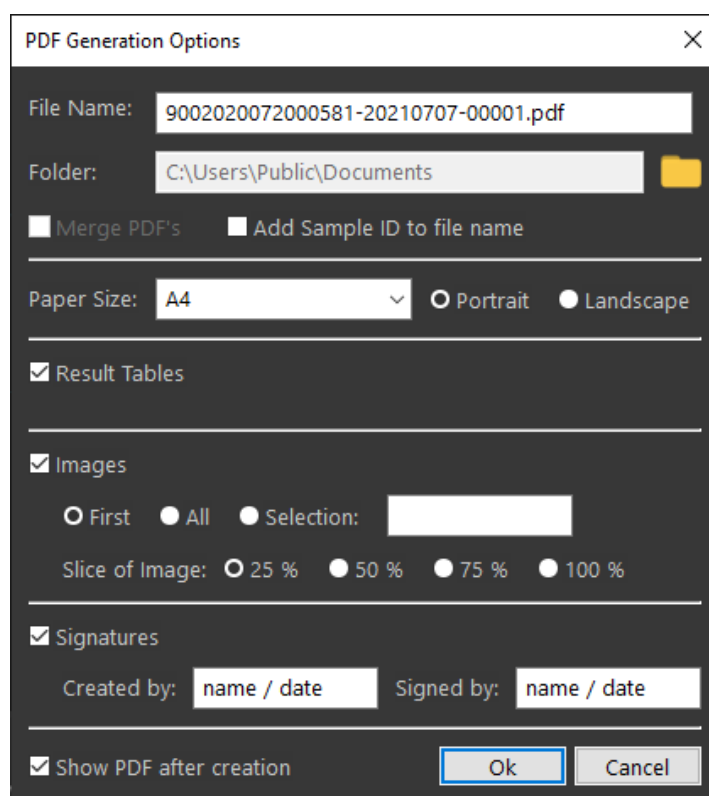


The screenshot shows the 'PDF Generation Options' dialog box. It has a title bar with a close button. The 'File Name' field contains '9002020072000581-20210707-00001.pdf'. The 'Folder' field shows 'C:\Users\Public\Documents' with a folder icon to its right. Below these are two checkboxes: 'Merge PDFs' (unchecked) and 'Add Sample ID to file name' (unchecked). The 'Paper Size' section has a dropdown set to 'A4' and two radio buttons for 'Portrait' (selected) and 'Landscape'. There are three checked checkboxes: 'Result Tables', 'Images', and 'Signatures'. Under 'Images', there are radio buttons for 'First' (selected), 'All', and 'Selection:' (with an empty text field). Below that are radio buttons for 'Slice of Image' with options '25 %' (selected), '50 %', '75 %', and '100 %'. Under 'Signatures', there are two text fields: 'Created by: name / date' and 'Signed by: name / date'. At the bottom, there is a checked checkbox 'Show PDF after creation' and two buttons: 'Ok' and 'Cancel'.

Software Options

Locate software options in the menu line of the [Main Window](#). Select *Tools* in the menu line and then *Options*. The Options window has three tabs: *General*, *Auto Export* and *Authorization*. In the *Auto Export* tab, the application can be setup to autogenerate CSV and/or PDF reports for different events. From the *Authorization* tab, NC-View™ can be configured to require a user to log in when launching the application. Furthermore, this tab can be used to activate the *Secure Mode* feature. For details see the NC-View™ NC-202™ Secure Mode Guide (doc. No. 991-2021).

Auto-export CSV files


1. To auto-export CSV files, select the *Auto Export* tab in the Options window.

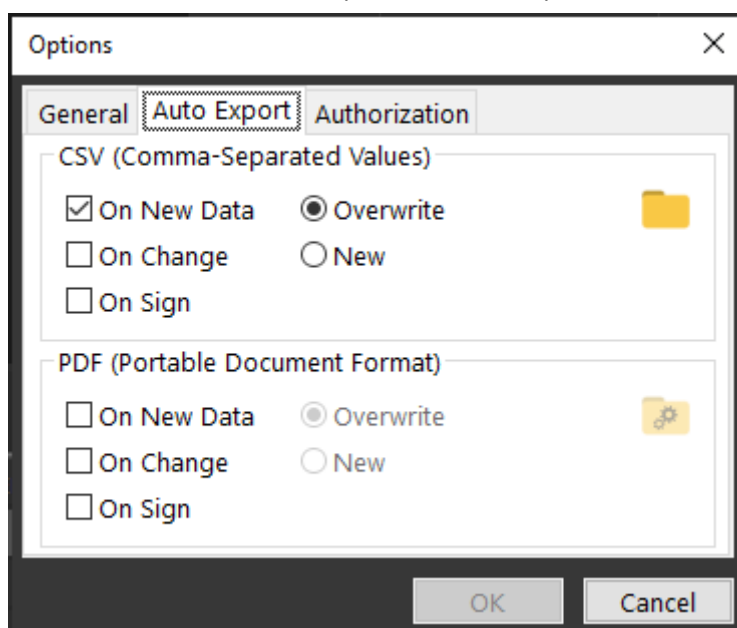

2. Select the events that trigger the auto-export of CSV files; select 'On New Data' to enable auto-export when new data is acquired; select 'On Change' to enable auto-export when data is changed (this includes changing the Sample ID, adding or deleting a tag, or adding a comment to a data set); select 'On Sign' to enable auto-export when data is signed.
3. Choose whether the CSV file should be overwritten with a new event or if a new CSV file should be created with every new event enabled.
4. Click the folder icon to select the destination folder.

Auto-export PDF reports

1. To auto-export PDF files, select the *Auto Export* tab in the Options window

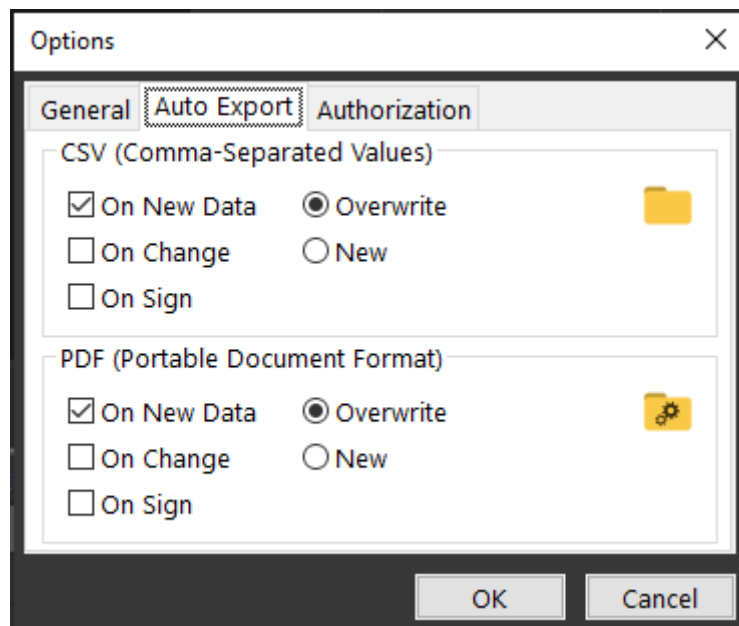

2. Select the events that trigger auto-export PDF reports; select 'On New Data' to enable auto-export when new data is acquired; select 'On Change' to enable auto-export when data is changed (this includes changing the Sample ID, adding or deleting a tag, or adding a comment to a data set); select 'On Sign' to enable auto-export when data is signed.
3. Select whether the PDF report should be overwritten with a new event, or if a new PDF report should be created with every new event enabled.
4. Click the folder/setup icon to open the PDF Generation Options and select destination folder, paper size, Result Tables, Images and Signatures.

License(s) and protocol limitations

NucleoCounter® NC-202™ instruments are sold with various licenses, e.g. the Secure Mode license, which allows acquisition of data in Secure Mode. Protocols available are restricted to the license(s) installed.

Certificate of Product Testing

When connected to an instrument, the Certificate of Product Testing for the instrument can be opened using the *About* box. This is opened by selecting *Help* in the menu line of the [Main Window](#) followed by *About*.

Installation Guide

Computer Specifications for Running NC-View™ Software:

Recommended minimum hardware and software requirements for running the NC-View™:

- Operating System: Windows 10 or 11. Prior to installing NC-View™, install the latest service packages and any critical updates to the operating system.
- Windows administrator credentials are required to install NC-View™.
- Processor: Intel Core i7 or i9 (9th or 10th generation); 2.5 GHz frequency (clock speed) or more.
- Hard drive: M2 SSD disk (500 GB) and at least 50 GB free disc space.
- RAM: 16 GB RAM or more.
- USB: USB 3.1 port (type A or type C).
- Display: 1920 × 1080 pixels screen resolution.

6.

NOTE: Computers with these specifications will perform a Count & Viability assay in 25-30 seconds. A lower CPU frequency will increase analysis time. For cleanroom applications requiring tablet computers, we recommend Microsoft Surface Pro Intel® Core™ i5 processor and 8 GB RAM, which will perform a Count & Viability assay in 30-37 seconds. NC-View™ installation will proceed regardless of computer specifications, but Windows 10 or 11 OS, USB 3.1, and administrator rights are a requirement.

Installing and Upgrading

If the NC-View™ application has previously been installed, the previous version will be removed during a new installation. Acquired data and saved protocols will not be deleted.

1. Ensure that the NC-202™ instrument is disconnected from the computer.
IMPORTANT: Log on with administrator rights for the installation session.
2. Open the location of the installation files in Windows File Explorer.
3. Double-click on the *Install NC-View X.X.X.X.exe* file to launch the installation program. The version number will vary depending on the software build, e.g. 2.1.0.21.
NOTE: Double-click on the .exe file, NOT the .bin file.
4. Follow the instructions until the installation is complete.
NOTE: During the installation, dialogs for installing third-party drivers will appear (e.g. Point Grey camera and other device drivers). Installing these drivers is required.

Uninstall Procedure

1. Right-click on *NC-View* in the Windows start menu and select *Uninstall*.
2. Follow the instructions until the uninstall process is completed.

NOTE: Uninstalling NC-View™ does not delete protocols and acquired data.

Importing protocols

1. In the menu line of the [Main Window](#), select *File* and then *Import Package*.
2. Locate the path where the protocol package is stored and select *OK* to import the protocol(s).

Maintenance and Backup

A CM file is approximately 25 MB in size. Consequently, acquiring many CM files will accumulate a large amount of data. We therefore recommend to backup CM files and free up disk space in good time to ensure trouble-free operation.

To export bulk amounts of CM files from a specific time period, open the [File Browser](#) and select the time period for which backup is required. Select all files by using the 'ctrl' and 'a' keys. Drag the selected files to an external hard disk or a network drive.

When backup is secured, the backed-up files can be deleted from the File Browser. Be aware that NC-View™ will add a comment to the CM files if they are exported or copied to show that the data integrity may have been compromised.

Troubleshooting

For support and troubleshooting regarding the NucleoCounter® NC-202™ and NC-View™ please contact ChemoMetec Support at support@chemometec.com

In case the solutions below do not solve the issue at hand, please create [Support Files](#) and contact ChemoMetec Support.

Issue	Solutions/options
Installation <i>The NC-View™ software or the NucleoCounter® NC-202™ instrument is not installed correctly.</i> <i>Server (ChemoMetec service) error.</i> <i>Camera not connected via USB3.</i>	<p>The user must be logged on with administrator rights during installation to enable correct installation.</p> <p>The installation can be blocked if an update process is running in the background, <i>e.g.</i> Windows update.</p> <p>If the installation of NC-View™ fails, ensure the Windows update is complete and try to run the installation process again using administrator rights login.</p> <p>NC-View™ requires a ChemoMetec server to run automatically after startup of the computer.</p> <p>Please restart the computer.</p> <p>A warning is displayed if the USB connection to the instrument initializes as a USB 2 connection. Please ensure the following:</p> <ol style="list-style-type: none">1) The USB cable supplied by ChemoMetec is used.2) A USB 3.1 port on the computer is used.3) The USB cable is securely screwed into the instrument.4) The USB cable is correctly inserted into the USB 3.1 port on the computer. Make sure that the USB cable is not bent in the USB 3.1 port.5) The cable is connected directly to the computer (<i>e.g.</i> do <u>NOT</u> use USB extender, adaptor or hub).6) If the computer is a desktop computer, use a USB 3.1 port on the back of the computer.7) Test a different USB 3.1 port on the computer to see if this resolves the issue.

<p>NC-View™ software</p> <p><i>Export of results data with the wrong decimal separator.</i></p>	<p>When copying result data from the file browser or file list, NC-View™ will place the data on the clipboard with the regional setting selected in the Windows OS.</p> <p>Be aware that if the region format has been set to <i>Match Windows display language</i>, changing the Windows decimal separator symbol will have no effect.</p> <p>Some applications, like spreadsheet software applications, also have their own options for setting the decimal separator symbol.</p>
<p>Via2-Cassette™</p> <p><i>No cassette is detected.</i></p> <p><i>Flow detection error.</i></p> <p><i>Cassette error.</i></p> <p><i>The cassette is stuck in the instrument.</i></p>	<p>This may be because either no Via2-Cassette™ is detected in the instrument or the Via2-Cassette™ is placed incorrectly. In the latter case, remove the Via2-Cassette™ and re-insert it.</p> <p>The sample loaded into the Via2-Cassette™ has failed to reach the counting chamber within the given time limits. Inserting an empty or an already used Via2-Cassette™ will trigger this error. If the sample volume is too small, it may also give this error.</p> <p>The inserted cassette is already used or is not clean. Take care not to touch the cassette windows while handling cassettes.</p> <p>Click <i>Ok</i> if any dialog boxes pop-up with this error. If the LED indicator on the instrument is green, click on the <i>Eject</i> icon. This will position the piston motor correctly.</p> <p>If this does not help or the LED indicator is red or orange, turn off the instrument via the main switch or socket, wait for ten seconds and then turn it back on. Check the USB connection between the instrument and the computer.</p>

