

# Chapter1 OverView

## 1.6 Measurement Samples

---

The TD-20's measurement sample is a sample tube of length 90 mm and outer diameter 1/4". The following types of sample tube are available. No sample tube is provided with the TDTS-2010. Obtain one separately.

223-52284-91	Glass sample tube: Tenax GR (130 mg)
223-52884-91	Glass sample tube: Tenax TA (100 mg) + Carboxen 1000 (50 mg)
223-57102-91	Glass sample tube: Tenax TA (130 mg)
223-57119	Glass sample tube (empty)
223-57474-91	Glass sample tube Carbopack(130 mg) + Carboxen 1000 (100 mg)

Special caps are attached to both ends of the sample tube and the tube is set in sample tray. Caps are not provided. Obtain them separately.

223-54617-91	Caps (20 per set)
--------------	-------------------

### **CAUTION**

Using a sample tube other than the ones specified may result in damage to the equipment.